



AN 382 Quantitative Damage Profiles By Ion Channeling

May 7, 2007 (Version 4.0)

Discussion

Ion channeling is used to study crystal damage, defect concentration, epitaxial crystal alignment, etc. When an ion beam is aligned along a major crystal axis or plane, ion-atom interaction probability is significantly reduced resulting in a large reduction of scattering events and deeper penetration of ions into the crystal. Simultaneously, redistribution of the ion flux in the channel occurs resulting in a phenomenon called flux-peaking. The ion flux is higher in the center of the channel than near the crystal axis or plane. Consequently, defects which are not located perfectly on the crystal lattice have a larger probability of scattering incoming ions.

In the backscattered spectrum from a channeled sample there are two components: the channeled ions scattered by defects in the crystal, and the de-channeled ions scattered by all atoms in the crystal. This method extracts the component of the channeled ions scattered from the defects and provides a quantitative damage profile in the crystal.

Figure 1 shows the channeling yield raw data plotted versus depth from a Si sample implanted with $1E14$ atoms/cm² of P. The extracted quantitative damage profile is shown in Figure 2. For comparison the simulated damage profile obtained from a TRIM simulation is shown in Figure 3. The TRIM damage profile shows the total damage (displacement) for all atoms in an amorphous Si sample. The sample shown in Figures 1 and 2 had P implanted in a channeling direction; the real damage profile is shallower than in the simulated amorphous sample.

Other Applications

- Quantitative impurity lattice locations (interstitial or substitutional) in crystal: Li to U by RBS(Rutherford Backscattering), PIXE(Particle Induced X-ray Emission) and NRA (Nuclear Reaction Analysis) with channeling
- Quality of Epitaxial Layer: Defect and Dislocation

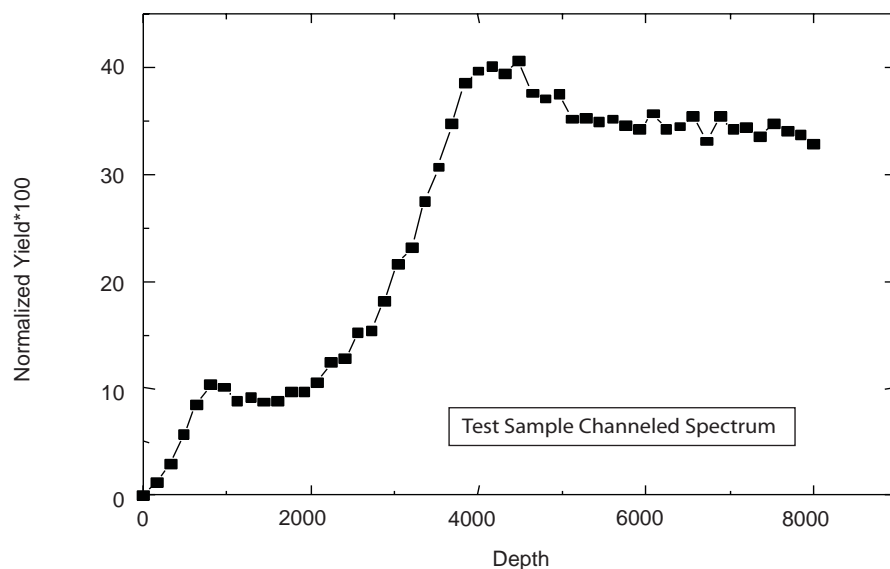


Figure 1. Channeling Yield Raw Data

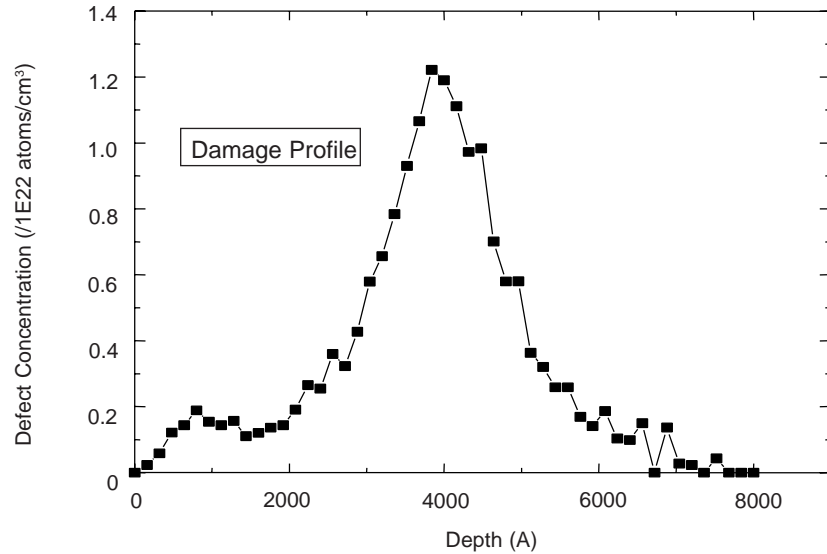


Figure 2. Quantitative Damage Profile

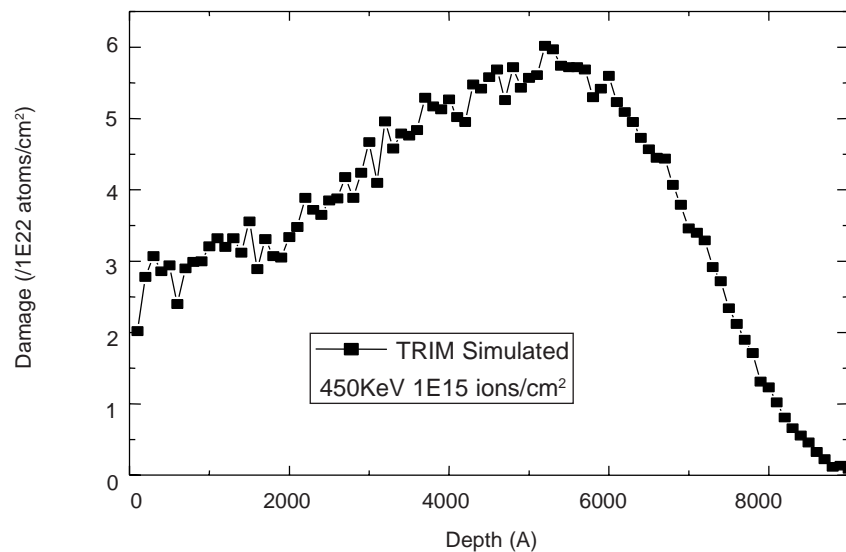


Figure 3. TRIM Simulated Damage Profile in Amorphous Si

United States Locations

Tempe, Arizona
+1 480 239 0602 info.az@eaglabs.com
+1 602 470 2655 fax

Sunnyvale, California
810 Kifer Road
+1 408 530 3500 info.ca@eaglabs.com
+1 408 530 3501 fax

1135 E Arques Avenue
+1 408 738 3033
+1 408 738 3035 fax

785 Lucerne Drive
+1 408 737 3892
+1 408 737 3916 fax

Peabody, Massachusetts
+1 978 278 9500 info.ma@eaglabs.com
+1 978 278 9501 fax

Chanhassen, Minnesota
+1 952 828 6411 info.mn@eaglabs.com
+1 952 828 6449 fax

East Windsor, New Jersey
+1 609 371 4800 info.nj@eaglabs.com
+1 609 371 5666 fax

Syracuse, New York
+1 315 431 9900 info.ny@eaglabs.com
+1 315 431 9800 fax

Raleigh, North Carolina
+1 919 829 7041 info.nc@eaglabs.com
+1 919 829 5518 fax

Round Rock, Texas
+1 512 671 9500 info.tx@eaglabs.com
+1 512 671 9501 fax

International Locations

Shanghai, China
+ 86 21 6879 6088 info.cn@eaglabs.com
+ 86 21 6879 9086 fax

Tournefeuille, France
+ 33 5 61 73 15 29 info.fr@eaglabs.com
+ 33 5 61 73 15 67 fax

Frankfurt, Germany
+ 49 (0) 693053213 info.de@eaglabs.com
+ 49 (0) 69307941 fax

Tokyo, Japan
+ 81 3 5396 0531 info.jp@eaglabs.com
+ 81 3 5396 1930 fax

HsinChu, Taiwan
+ 886 3 5632303 info.tw@eaglabs.com
+ 886 3 5632306 fax

Uxbridge, United Kingdom
+ 44 (0) 1895 811194 info.uk@eaglabs.com
+ 44 (0) 1895 810350 fax